

USP <800> Surface Wipe Sampling Test

Let Us Help You Verify Workplace Safety in 4 Easy Steps

1 Buy Texwipe Swabs Part No TX714K

Texwipe swabs enhance absorbency, increase lifting, and capture small particles.

**Available through
Letco Med**

Product ID: 696632

Call (800) 239-5288 or visit
letcomed.com to order today

2 Identify Swab Areas

According to USP <800>, swab sampling should be performed routinely.

Sample Collection Areas:

- Inside of C-PEC
- Pass-through chambers
- Staging and work surfaces
- Floors and dispensing areas
- Outside HD buffer room or C-SCA
- Patient administration

3 Collect Sample

- Wet swab and shake off excess water
- Swab sampling area
- Place swab in vial
- Cut swab handle at perforation
- Cap Vial
- Label vial with sample collection area
- Change gloves and repeat process for other sampling areas

4 Submit Sample to ARL

Visit www.arlok.com/HDs for list of available drugs. Call ARL for drugs not listed.

Sample Shipment:

- Package vials for shipping and include cool packs
- Write USP <800> testing on sample submission form and list drugs to be tested
- Results delivered in 10 business days

\$200 for 1st HD + \$50 for each additional HD per swab sample

CONTACT ARL TODAY

Email: info@arlok.com

Website: www.arlok.com

Toll Free: (800) 393-1595